Pressure gauge for sanitary applications
Models PG43SA-S, PG43SA-C

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Prior to starting any work, read the operating instructions!
 Keep for later use!

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Contents

1. General information	4
2. Design and function	5
3. Safety	6
4. Transport, packaging and storage	7
5. Commissioning, operation	9
6. Faults	12
7. Maintenance and cleaning	12
8. Dismounting and disposal	13
9. Specifications	14

Declarations of conformity can be found online at www.wika.com.

1. General information

- The pressure gauge described in the operating instructions has been designed and manufactured using state-of-the-art technology. All components are subject to stringent quality and environmental criteria during production. Our management systems are certified to ISO 9001 and ISO 14001.
- These operating instructions contain important information on handling the instrument. Working safely requires that all safety instructions and work instructions are observed.
- Observe the relevant local accident prevention regulations and general safety regulations for the instrument's range of use.
- The operating instructions are part of the product and must be kept in the immediate vicinity of the instrument and readily accessible to skilled personnel at any time. Pass the operating instructions on to the next operator or owner of the instrument.
- Skilled personnel must have carefully read and understood the operating instructions prior to beginning any work.
- The general terms and conditions contained in the sales documentation shall apply.

Notes per pressure equipment directive

These instruments are defined as "pressure accessories" per pressure equipment directive chapter 1, article 2.5.

Applied standards

EN 837-2 Selection and installation recommendations for pressure gauges

EN 837-3 Diaphragm and capsule pressure gauges, dimensions, metrology, requirements and testing

2. Design and function

- Subject to technical modifications.
- Further information:
 - Internet address: www.wika.de / www.wika.com
 - Corresponding data sheets: PM 04.15, PM 04.16
 - Application consultant: Tel.: +49 9372 132-0
Fax: +49 9372 132-406
info@wika.de

EN

2. Design and function

2.1 Description

The model PG43SA-S and PG43SA-C pressure gauges are designed especially for the requirements of the food and beverage, pharmaceutical and biotechnology industries.

The pressure gauge is particularly suitable for the special conditions of CIP/SIP cleaning processes, such as chemical stability towards cleaning liquids and high temperatures. The flush metal diaphragm element is directly welded to the process connection. This guarantees a crevice-free connection of process connection and diaphragm element; additional internal sealings are not required.

2.2 Scope of delivery

Cross-check scope of delivery with delivery note.

3. Safety

3.1 Explanation of symbols

WARNING!

... indicates a potentially dangerous situation that can result in serious injury or death, if not avoided.

Information

... points out useful tips, recommendations and information for efficient and trouble-free operation.

3.2 Intended use

The model PG43SA-S and PG43SA-C pressure gauges are used for measuring pressure particularly in sanitary applications.

The instrument has been designed and built solely for the intended use described here, and may only be used accordingly.

The manufacturer shall not be liable for claims of any type based on operation contrary to the intended use.

3.3 Improper use

WARNING!

Injuries through improper use

Improper use of the instrument can lead to hazardous situations and injuries.

- ▶ Refrain from unauthorised modifications to the instrument.

Any use beyond or different to the intended use is considered as improper use.

3. Safety

EN

3.4 Compliance with the conformity in accordance with 3-A

For a 3-A compliant connection the following sealings have to be used:

- For milk thread fittings per DIN 11851, suitable profile sealings have to be used (e.g. SKS Komponenten BV or Kieselmann GmbH).
- For fittings per IDF sealings with support ring per ISO 2853 have to be used.

Note: Connections per SMS, APV RJT and NEUMO Connect S are not 3-A compliant.

3.5 Compliance with EHEDG conformity

For an EHEDG conform connection, sealings in accordance with the current EHEDG policydocument must be used.

Sealings for connections per ISO 2852, DIN 32676 and BS 4825 part 3 are, e.g., manufactured by Combifit International B.V. A manufacturer of sealings for connections per DIN 11851 is, e.g., Kieselmann GmbH. A manufacturer of VARIVENT® sealings is, e.g., GEA Tuchenhagen GmbH. A manufacturer of NEUMO BioConnect® sealings is, e.g., Neumo GmbH & Co. KG.

3.6 Personnel qualification

WARNING!

Risk of injury should qualification be insufficient
Improper handling can result in considerable injury and damage to equipment.

- ▶ The activities described in these operating instructions may only be carried out by skilled personnel who have the qualifications described below.

Skilled personnel

Skilled personnel, authorised by the operator, are understood to be personnel who, based on their technical training, knowledge of measurement and control technology and on their experience and knowledge of country-specific regulations, current standards and directives, are capable of carrying out the work described and independently recognising potential hazards.

4. Transport, packaging and storage

4.1 Transport

Check the measuring instrument for any damage that may have been caused by transport.

Obvious damage must be reported immediately.

CAUTION!

Damage through improper transport

With improper transport, a high level of damage to property can occur.

- ▶ When unloading packed goods upon delivery as well as during internal transport, proceed carefully and observe the symbols on the packaging.
- ▶ With internal transport, observe the instructions in chapter 4.2 “Packaging and storage”.

4.2 Packaging and storage

Do not remove packaging until just before mounting.

Keep the packaging, especially the diaphragm protection (if available). This will provide optimum protection during transport (e.g. cleaning, change in installation site, sending for repair).

CAUTION!

Damage to the diaphragm during unpacking

With improper unpacking, a high level of damage to property can occur.

Avoid touching the diaphragm after removing the diaphragm protection. Impermissible mechanical loading (e.g. pressing the diaphragm) can result in measuring deviations or even in the complete failure of the measuring instrument.

The PG43SA-C is optionally packed under sterile conditions, and on the perforation on the rear of the sterile packaging there is a brown-coloured process indicator. By checking this brown colouration you can find out whether the gauge has been packed under sterile conditions.

Store instruments that are packed under sterile conditions in the instrument package (no condensation allowed, storing possible for 6 months).

Permissible conditions at the place of storage

Storage temperature: -30 ... +70 °C

Protect the instrument from dust.

5. Commissioning, operation

5.1 Mounting preparation

- To protect the pressure gauge from mechanical damage keep it in the outer carton until installation.
- The gauge should only be removed from the sterile packaging shortly before mounting at the measuring point.

5. Commissioning, operation

EN

- With CIP (cleaning in place) we recommend to clean the measuring instrument again after mounting, using the process established at the operator's site.
- With COP (cleaning out of place) we recommend to clean the measuring instrument again before mounting, using the process established at the operator's site.
- On removal from the packaging, and during mounting, particular care must be taken to avoid damage and any mechanical deformation to the diaphragm.
- Model PG43SA-S only: Check the zero point. In case of a zero offset, carry out a zero point setting. See chapter 5.7 "External zero point setting".

5.2 Installation

- Do not damage the diaphragm; scratches on the diaphragm (e.g. from sharp-edged objects) are the main causes of corrosion.
- For sealing, choose appropriate seals.
- For flange mounting, only use sealings with a sufficiently large inner diameter and centre them; contact with the diaphragm leads to measuring errors.
- When using soft or PTFE sealings, observe the instructions of the sealing manufacturer, particularly with regard to tightening torque and load cycles.
- For installation, in accordance with the fitting and flange standards the appropriate fastenings, such as screws and nuts, must be used.
- The vent hole at the underside of the case must not be closed (only applies to model PG43SA-S).

5. Commissioning, operation

- Nominal position per EN 837-3 / 9.6.6 figure 7: 90° (⊥) (only applies to model PG43SA-S)

5.3 Mounting instructions for instruments with EHEDG and 3-A

Observe the following instructions, especially for EHEDG certified and 3-A conform instruments.

- To maintain the EHEDG certification, one of the EHEDG-recommended process connections must be used. These are marked with the logo in the data sheet.
- To maintain the conformity to the 3-A standard, a 3-A conform process connection must be used. These are marked with the logo in the data sheet.
- Mount the instrument with minimal dead space and able to be cleaned easily.
- The mounting position of the instrument, welding socket and instrumentation T-piece should be designed to be self-draining.
- The mounting position must not form a draining point or cause a basin to be formed.
- With the process connection via an instrumentation T-piece, the branch L of the T-piece must not be longer than the diameter D of the T-piece ($L \leq D$).

Special instructions for in-line versions

- In-line access units (e.g. NEUMO BioControl®) and instrumentation T-pieces in horizontal pipes should be installed with a slight pipe inclination for self-draining.
- With in-line access units (e.g. NEUMO BioControl®), use the original seals from the manufacturer and observe their installation instructions.

5.8 Autoclaving

The following applies only to model PG43SA-S with option “autoclavable version”, recognizable on the dial printing as “AUTOCLAVABLE”.

- Observe the permissible conditions of autoclaving, see chapter 9.3 “Permissible temperature”
- During the autoclaving process, no weight-loading must be exerted via the case and the instrument must not lie on its window.
- After the first autoclaving cycles, the window slightly arches outward. This has no effect on the function of the instrument.
- Depending on the autoclaving process, whitish residues can build up on the window and pointer. This has no effect on the function of the instrument.
- In case of a zero offset, carry out a zero point setting. See chapter 5.7 “External zero point setting”.

6. Faults

Following any failure or shut-down of the plant all instruments should be checked and, if necessary, replaced before recommissioning the plant.

7. Maintenance and cleaning

7.1 Maintenance

- The instruments are maintenance-free.
- The indicator should be checked once or twice every year. For this the instrument must be disconnected from the process to check with a pressure testing device.

- Repairs should only be carried out by the manufacturer or, following prior consultation with the manufacturer, by correspondingly qualified skilled personnel.

7.2 Cleaning

Suitable for 'in place' SIP and CIP cleaning processes, if appropriate media are used and the permissible specifications are complied with.

When cleaning from outside ("wash down"), observe the permissible temperature and ingress protection.

For sterilisation in an autoclave, observe the instructions in chapter 5.8 "Autoclaving".

Vent hole with model PG43SA-S

In order to avoid any change in pressure within the case, the model PG43SA-S features a vent hole at the case bottom on the rear side. In case of condensation formation within the case or ingress of low water quantities from outside, this can be drained through the vent hole.

8. Dismounting and disposal

WARNING!

Physical injuries and damage to property and the environment through residual media

Residual media in the instrument can result in a risk to persons, the environment and equipment.

- ▶ Wash or clean the dismantled instrument, in order to protect persons and the environment from exposure to residual media.

8. Dismounting and disposal / 9. Specifications

EN

8.1 Dismounting

Only disconnect the measuring instrument once the system has been depressurised!

8.2 Disposal

Incorrect disposal can put the environment at risk.

Dispose of instrument components and packaging materials in an environmentally compatible way and in accordance with the country-specific waste disposal regulations.

9. Specifications

9.1 Pressure limitation

Model	Steady	Fluctuating	
		Scale range < 4 bar	Scale range \geq 4 bar
PG43SA-S	Full scale value	Full scale value	2/3 x full scale value
PG43SA-C	3/4 x full scale value	2/3 x full scale value	

9.2 Overload safety

Model	Overload safety ¹⁾
PG43SA-S	2 x full scale value, max. 40 bar or max. pressure rating (PN) of the process connection
PG43SA-C	5 x full scale value, max. 40 bar or max. pressure rating (PN) of the process connection

1) see PN in the tables of data sheets PM 04.15, PM 04.16

9.3 Vacuum safety (only model PG43SA-S)

All scale ranges are suitable for the vacuum typical of CIP or SIP cleaning. Positive scale ranges < 6 bar (< 100 psi) are optionally available in a vacuum-tight version (to -1 bar).

9. Specifications

EN

9.4 Permissible temperature

Place, Situation	Permissible temperature
Ambient	-20 ... +60 °C
Medium	≤ 150 °C
CIP and SIP	150 °C continuously for wetted parts
Autoclaving ¹⁾	≤ 134 °C, 20 minutes
Storage, transport	-30 ... +70 °C

1) only for model PG43SA-S with option "autoclavable version"

9.5 Temperature effect

Model	Deviation from the reference temperature (+20 °C) at the measuring system
PG43SA-S	≤ ±0.5 %/10 K of full scale value
PG43SA-C	≤ ±0.8 %/10 K of full scale value

9.6 Ingress protection per IEC/EN 60529

Model PG43SA-S: IP54

Model PG43SA-C: IP66

For further specifications see WIKA data sheets PM 04.15, PM 04.16 and the order documentation.

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All rights reserved. / Alle Rechte vorbehalten.
WIKA® is a registered trademark in various countries.